Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The instrument is not suitable for the use with wood, stone and metal.
- ▶ Do not use the instrument on live humans or animals.
- ▶ Wear the appropriate protective clothing.

Layout and function

- ▶ The instrument is made of stainless steel 1.4034 (AISI 420).
- ▶ The instrument is suitable to take samples of semi-solid materials in the food and beverage industry using hand-held drills or cordless screwdrivers.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is maintenance-free and suitable for

steam sterilisation (up to 121°C) for sterilization by flame treatment.

- ▶ If necessary, the cutting edge can be resharpened.

Parts in contact with medium

- ▶ Stainless steel 1.4034 (AISI 420).

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Mount the drill head ① on a hand-held drill or cordless screwdriver.

NOTE!

Verify that the borer is attached correctly before use.

Safe sampling

WARNING!

Bodily injury due to accidental contact with electrical power. Danger of electric shock.

- ▶ Keep the cable of the drill out of the cutting area.

WARNING!

Sharp parts on the instrument. Risk of injury.

- ▶ Keep hands well away from the cutting area.

WARNING!

Bodily injury due to drilling through the sampling material without adequate securing.

- ▶ Use a suitable base (e.g. cutting board made of plastic or wood).
- ▶ Use a suitable locking mechanism or secure the sampling material against twisting and shifting.
- ▶ Only place the instrument vertically onto the sampling material.

WARNING!

Bodily injury due to entanglement of body parts, hairs or clothes.

- ▶ Keep hands well away from the cutting area.
- ▶ Use a hairnet if necessary.
- ▶ Ensure to wear tight-fitting clothes.

1. Place the drill head ① onto the sampling material.
2. Start boring slowly until a slight trace is incised in the sampling material.
3. Pull the drill head ① out of the sampling material.
4. Remove the sample carefully from the sampling material.
5. Clean the instrument.

CAUTION!

Sharp parts on the instrument. Risk of injury.

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